

VeraMag™

REF 500020 and 500021

Email: info@veravas.com



Intended Use

The Veravas VeraMag™ is a magnetic particle separation device that uses neodymium magnets to separate different biomolecules for further analysis.

SUMMARY AND EXPLANATION

VeraMag™ supports a complete range of magnetic separation applications including VeraPrep Biotin™ (Part No. 500014 and 500015), RNA and DNA isolation, and biomolecules purification. It consists of an injection molded plastic housing which incorporates a neodymium magnet.

It is designed to accommodate micro tube 2ml with cap (SARSTEDT Order Number 72.694), standard 1.5 ml Conical tubes, and standard 2.0 mL HPLC tubes.

REAGENTS AND MATERIALS PROVIDED

CONTENT		
DEVICE	Magnetic particle separation device	
REF	500020	500021
DEVICE	1	2

MATERIALS REQUIRED BUT NOT PROVIDED

1. Pipetting device(s) capable of delivering 50 µL up to 1000 µL
2. Disposable pipette tips
3. Micro tube 2ml with cap (SARSTEDT Order Number 72.694), standard 1.5 ml Conical tubes, or standard 2.0 mL HPLC tubes
4. Personal protective equipment

STORAGE AND STABILITY

Upon receipt, store in the box at 5°- 25°C.

Warnings and Handling

- EXPORT**
1. Veravas VeraMag products should be handled with care. Avoid direct contact with the magnet material. Persons with pacemakers/implants should avoid direct contact. Keep loose ferrous material away and do not attempt to disassemble.

VERAVAS

transforming laboratory medicine

2. Keep all magnetic media, watches, and sensitive electronic devices away from the units.
3. Credit cards, tape and disks can be erased in the presence of a magnetic field. Bodily harm (pinching of hands and skin) can result if magnets are not handled correctly. Maintain distance between two or more units.
4. The product should be stored in a cool, dry environment. Clean using a damp, lint-free cloth. Sanitize using a mild detergent, (10% Bleach Solution or IPA).
5. Do not spray solution directly onto unit. Spray on cloth and wipe clean. Sani-Cloth Wipes may also be used.
6. Do not autoclave.



Qarad EC-REP BV
Pas 257, 2440 Geel
BELGIUM



Veravas Inc
128 Columbus Street, # 825,
Charleston, SC, 29403 USA



DEVICE

VeraMag™ magnetic particle separation device

EXPORT

For Export Only